

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Illumination pulsing on the monitor	Bulletin No:	SB0914
	Date Issued:	6/25/98
	Priority:	Medium

Product: LSX-200

Serial Numbers: ALL
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

The LSX-200 monitor exhibits illumination pulsing that may indicate it is out of sync.

IMPLEMENTATION

If the monitor seems to be flashing the problem may be the potentiometer on the dual lamp supply board in the LSX-200. The potentiometer may need to be adjusted. If the LSX-200 is equipped with a GEO option verify by looking through the binoculars that the problem is indeed with the monitor being out of sync. If the flashing is only occurring on the monitor remove the panel on the right side of the LSX-200 and locate the lamp supply board. Adjust the potentiometer until the monitor stops flashing.

PARTS

Part Number	Description
-------------	-------------

AVAILABLE DOCUMENTATION