

# CETAC TECHNOLOGIES

## SERVICE BULLETIN

<b>Subject:</b> Main cabinet does not power-up	<b>Bulletin No:</b>	SB0915
	<b>Date Issued:</b>	6/26/98
	<b>Priority:</b>	Medium

**Product:** LSX-200

**Serial Numbers:** ALL

*Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)*

**Service Location:**  Customer Site  CETAC Technologies

**Responsibility:**  Customer  Trained CETAC Representative

**Replacement Parts:**  Customer  CETAC Technologies

### DESCRIPTION

The LSX-200 main cabinet will not power-up.

---

---

---

---

### IMPLEMENTATION

If the main cabinet does not power-up or initialize the problem may be a blown fuse. If the fuses ohm out ok the supply voltage at the power supply needs to be verified. If power exists at the power supply and the main cabinet does not power-up, then the problem may lie within the power supply. Unplug the power out of the power supply so the voltage can be checked at the power supply output. The power out leads on the power supply should be supplying 24V DC. If not, then the power supply is the problem..

---

---

### PARTS

Part Number	Description
-------------	-------------

### AVAILABLE DOCUMENTATION