

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Autosampler X-Axis Sensor Board Replacement	Bulletin No:	SB1001
	Date Issued:	9/21/2004
	Priority:	Medium

Product: ASX-510HS, ASX-520

Serial Numbers: 470976 on X-axis board
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Replacement of X-axis sensor board to upgrade 4-pin connector to 6-pin connector.

IMPLEMENTATION

Replace X-axis sensor by removing the two screws and putting the new sensor in its place.

The Y-axis sensor also needs to be replaced. See available documentation for instructions.

Connect the Y-axis sensor to the X-axis sensor.

PARTS

Part Number	Description
SP6107	Y-axis sensor board
SP6106	X-axis sensor board

AVAILABLE DOCUMENTATION

SVC119.pdf Guide to Replacing the Y-Axis Sensor Board on the ASX-520 & ASX-520HS Autosamplers