

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Autosampler Power Connector Defective	Bulletin No:	SB1003
	Date Issued:	9/21/04
	Priority:	Low

Product: ASX-260, ASX-410, ASX-510HS, ASX-520
Serial Numbers: Prior to 100401A260, 100401A410, 100401A520
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Autosampler power connector has an inner connector that is weak that can break in shipping or when pressure is applied to it.

IMPLEMENTATION

Replace connector by removing screws and cutting the wires. Then solder wires to the new connector and putting the screws back on.

PARTS

Part Number	Description
SP6104	Switch/power input kit

AVAILABLE DOCUMENTATION