

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Elimination of “splashing” in high-speed autosamplers	Bulletin No: SB 1009
	Date Issued: 11/01/05
	Priority: Medium

Product: High-Speed Autosamplers

Serial Numbers: Before 090501A520HS

Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Quick ramp up of autosampler causes “splashing” of samples.

IMPLEMENTATION

Replace the main board with correct rabbit board attached using the instructions in the available documentation.

PARTS

Part Number	Description
SP7091	Main board with rabbit board(various versions) and installation guide.

AVAILABLE DOCUMENTATION